



EV231:	300 V,	1000 mA,	150 W
EV202:	300 V,	2000 mA,	300 W
EV245:	400 V,	500 mA,	50 W
EV265:	600 V,	500 mA,	150 W
EV261:	600 V,	1000 mA,	300 W
EV215:	1200 V,	500 mA,	300 W
EV233:	3000 V,	300 mA,	300 W
EV262:	6000 V,	150 mA,	300 W

Manual or Method programming
Computer control using RS232
Stores up to 3600 output values

Manual programming

The manual mode allows to set voltage, current, power and time for a routine electrophoresis run. Parameters can be changed temporarily without interrupting the run.

Method programming

Up to 9 different programs, each with 9 steps, of frequently used parameters can be stored in the non-volatile memory for future recall. Reads voltage in 1 V steps, current in 1 mA steps and power in 1 W steps. Each step is able to recall a next one, providing a flexible multiple step function for special techniques. Parameters of the running step can be changed temporarily without interrupting the run.

Voltage ramp

The method mode also allows to program a linear voltage gradient for any step provided the limiting current or power is not attained.

Timer

Timer or volt-hour controlled operation is a useful standard feature on all models. The microcomputer will automatically terminate the run and sound an alarm when the count down of the selected value is achieved.

Automatic cross-over

Each model has constant voltage, constant current, constant power capabilities with automatic cross-over and shows which parameter is kept constant.

Password protection

Prevents any unauthorised modification of the instrument's settings allowing the user only to run or stop an electrophoresis.

Automatic recovery after power failure

After a mains power failure the instrument will automatically continue the run for the remaining time.

Display

The interactive LCD screen provides step by step instructions in the language of your choice (English, Dutch, French, German).

Data-logging

Stores up to 3600 output values (voltage, current and power) including program number and step.

Data Transfer

Free data acquisition software for PC can be downloaded from our website. It allows to visualize and examine the stored run details via RS232.

Remote control

All power supplies can be controlled by a computer using special commands.

Safety precautions

The user is protected from potential shock hazard since the AC line is automatically disconnected from the high voltage transformer when a ground leakage path is detected.

The instrument is fully protected against any overload condition including accidental short circuit of the output.

The high voltage cannot suddenly appear at the outputs. It will always increase smoothly up to one of the pre-set limits is reached.

Galvanically isolated RS232 input/output connection prevents ground loop interferences when connected to a computer.

Warranty

Three year warranty.

Comprises: manual and mains lead.

SPECIFICATIONS	EV231	EV202	EV245	EV265
VOLTAGE	0...300 V	0...300 V	0...400 V	0...600 V
CURRENT	0...1000 mA	0...2000 mA	0...500 mA	0...500 mA
POWER	0...150 W	0...300 W	0...50 W	0...150 W
PARAMETER RANGE	1...100% of full scale	1...100% of full scale	1...100% of full scale	1...100% of full scale
TIMER	0...99:59 h	0...99:59 h	0...99:59 h	0...99:59 h
VOLT-HOURS	0...99.99 kWh	0...99.99 kWh	0...99.99 kWh	0...99.99 kWh
DISPLAY	LCD, 2x16 characters	LCD, 2x16 characters	LCD, 2x16 characters	LCD, 2x16 characters
RESOLUTION	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W
PROGRAMS	9x9 set of parameters	9x9 set of parameters	9x9 set of parameters	9x9 set of parameters
OUTPUTS	4 in parallel, 4 mm sockets	4 in parallel, 4 mm sockets	3 in parallel, 4 mm sockets	4 in parallel, 4 mm sockets
MINIMUM LOAD RESISTANCE	10 Ω	5 Ω	30 Ω	30 Ω
NO LOAD DETECTION	on/off, programmable	on/off, programmable	on/off, programmable	on/off, programmable
GROUND LEAKAGE DETECTION	✓	✓	✓	✓
OVERLOAD DETECTION	✓	✓	✓	✓
COMPUTER CONTROL	✓	✓	✓	✓
PASSWORD PROTECTION	✓	✓	✓	✓
DATA-LOGGING	3600 values	3600 values	3600 values	3600 values
INTERVAL	1...60 seconds	1...60 seconds	1...60 seconds	1...60 seconds
RS232 INTERFACE	9600 b/s	9600 b/s	9600 b/s	9600 b/s
AMBIENT TEMPERATURE	0...40°C	0...40°C	0...40°C	0...40°C
RELATIVE HUMIDITY	0...95%, non condensing	0...95%, non condensing	0...95%, non condensing	0...95%, non condensing
POWER REQUIREMENTS	210...250 VAC, 50/60 Hz, 200 W	210...250 VAC, 50/60 Hz, 360 W	210...250 VAC, 50/60 Hz, 75 W	210...250 VAC, 50/60 Hz, 200 W
DIMENSIONS (WxDxH)	31x26x15 cm	31x26x15 cm	24x20x13 cm	31x26x15 cm
WEIGHT	6 kg	10 kg	3 kg	6 kg

SPECIFICATIONS	EV261	EV215	EV233	EV262
VOLTAGE	0...600 V	0...1200 V	0...3000 V	0...6000 V
CURRENT	0...1000 mA	0...500 mA	0...300 mA	0...150 mA
POWER	0...300 W	0...300 W	0...300 W	0...300 W
PARAMETER RANGE	1...100% of full scale	1...100% of full scale	1...100% of full scale	1...100% of full scale
TIMER	0...99:59 h	0...99:59 h	0...99:59 h	0...99:59 h
VOLT-HOURS	0...99.99 kWh	0...99.99 kWh	0...99.99 kWh	0...99.99 kWh
DISPLAY	LCD, 2x16 characters	LCD, 2x16 characters	LCD, 2x16 characters	LCD, 2x16 characters
RESOLUTION	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W	1 V, 1 mA, 1 W
PROGRAMS	9x9 set of parameters	9x9 set of parameters	9x9 set of parameters	9x9 set of parameters
OUTPUTS	4 in parallel, 4 mm sockets	4 in parallel, 4 mm sockets	4 in parallel, 4 mm sockets	4 in parallel, 2 mm sockets
MINIMUM LOAD RESISTANCE	15 Ω	70 Ω	600 Ω	1200 Ω
NO LOAD DETECTION	on/off, programmable	on/off, programmable	on/off, programmable	on/off, programmable
GROUND LEAKAGE DETECTION	✓	✓	✓	✓
OVERLOAD DETECTION	✓	✓	✓	✓
COMPUTER CONTROL	✓	✓	✓	✓
PASSWORD PROTECTION	✓	✓	✓	✓
DATA-LOGGING	3600 values	3600 values	3600 values	3600 values
INTERVAL	1...60 seconds	1...60 seconds	1...60 seconds	1...60 seconds
RS232 INTERFACE	9600 b/s	9600 b/s	9600 b/s	9600 b/s
AMBIENT TEMPERATURE	0...40°C	0...40°C	0...40°C	0...40°C
RELATIVE HUMIDITY	0...95%, non condensing	0...95%, non condensing	0...95%, non condensing	0...95%, non condensing
POWER REQUIREMENTS	210...250 VAC, 50/60 Hz, 360 W	210...250 VAC, 50/60 Hz, 360 W	210...250 VAC, 50/60 Hz, 360 W	210...250 VAC, 50/60 Hz, 360 W
DIMENSIONS (WxDxH)	31x26x15 cm	31x26x15 cm	31x26x15 cm	31x26x15 cm
WEIGHT	10 kg	10 kg	10 kg	10 kg

CODE	DESCRIPTION
EV231	Power supply, 300 V, 1000 mA, 150 W
EV202	Power supply, 300 V, 2000 mA, 300 W
EV245	Power supply, 400 V, 500 mA, 50 W
EV265	Power supply, 600 V, 500 mA, 150 W
EV261	Power supply, 600 V, 1000 mA, 300 W
EV215	Power supply, 1200 V, 500 mA, 300 W
EV233	Power supply, 3000 V, 300 mA, 300 W
EV262	Power supply, 6000 V, 150 mA, 300 W
➔ Supplied with a european mains cord + USB cable	
(Add a US-sign for US plug 120 VAC versions, e.g.: EV245-US)	
(Add a UK-sign for UK plug versions, e.g.: EV245-UK)	
(Add a CH-sign for Swiss plug versions, e.g.: EV245-CH)	

Accessories

CODE	DESCRIPTION
E200	Pair of adaptors, 4 mm plug to 2 mm socket
E201	Pair of cables M/F, 4+4 mm
E204	Pair of adaptors, 2 mm plug to 4 mm socket
AK2315	RS232 comuter cable

Free data acquisition software for PC can be downloaded from www.consort.be

